

# Reduce Labor Costs and Increase Productivity with the CERTUS 3-Step Workflow

## Sample, Test, Result.

No specialized training, certification, or experience necessary.

Minimal prep steps means reduced hands-on time per sample.

Automation reduces repetitive manual tasks.


Now anyone in your facility can perform environmental pathogen testing. Simply collect your sample and insert into the Detection Unit following step-by-step directions on your Control Pad.



### 1 Sample



Collect sample using Bio-Lock™ Sampling Swab.

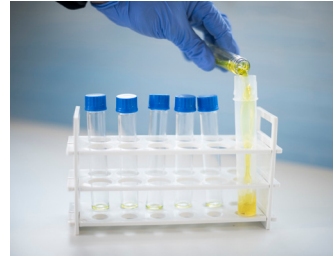

Pour pre-measured selective growth media into Detection Tube.



Insert Bio-Lock Sampling Swab into Detection Tube.

**BIO-LOCK™** Once securely closed, Bio-Lock Sampling Swabs form a permanent seal that can't reopen, thus securing the facility from accidental contamination.

### 2 Test



Initiate testing by clicking on "New Test" on Control Pad.



Scan the barcode on the Detection Kit box.



Insert Detection Tube into available slot.

### 3 Result



Real-time detection begins immediately with results in as little as 8 hours.



Test results on Detection Unit, Control Pad, and via email/SMS alerts.



All testing information is automatically logged and stored for easy traceability.